

*Esco PCR Cabinet, Model PCR-4A.
Shown with optional support stand.*

Polymerase Chain Reaction Cabinets *The Proven Solution for Contaminant-Free PCR*





Esco PCR Cabinet, Model PCR-4A_, shown with optional support stand.



Main Features

- Greater protection against contamination from the ambient environment and cross-contamination within the main chamber.
- High quality polyester pre-filter and main HEPA filter with a typical efficiency of $\geq 99.99\%$ at 0.3 microns.
- Built-in UV lamp with timer to facilitate decontamination between PCR cycles.
- Sentinel™ Microprocessor controller supervises all functions - 0.9 m (3') and 1.2 m (4') models only.
- Esco antimicrobial coating on all painted surfaces.
- Available in 0.6, 0.9 and 1.2 meter models (2', 3' and 4').



Introduction

What is PCR

Polymerase Chain Reaction (PCR) is a process where millions of copies of DNA are amplified from a single copy or low copy number template. This reaction is fundamental to almost all applications requiring a high copy number of starting material and is used in all laboratories

Why PCR Cabinet

Because of the high copy number generated and the nature of replication during PCR, it is essential to prevent possible contamination of the PCR reaction.

The ideal PCR laboratory should consist of three areas, each isolated from the other. Reagents should be prepared in the reagent preparation area and transferred to the sample preparation area through a pass box or inside closed containers. After preparation of the final reaction mix, the tubes should be transferred to the amplification area, again through a pass box or in a closed container. The PCR amplification and results analysis take place in this area.




In practice, these procedures may take place in the same room. Under this circumstances, PCR Cabinets are used for reagent preparation and sample preparation to minimize contamination. Where there remains a biological risk from the sample to the user and/or environment, Class II biological safety cabinets must be used for sample preparation.

How Esco PCR Cabinets Prevent Contamination

UV Decontamination Technology


- Powerful, uniformly distributed, UV-C.
- Proven effect on DNA contamination.
- Shelf enables placement of items closer to the UV source, increasing decontamination efficacy.
- UV-C is ozone-free.
- UV hour meter monitors bulb life, simplifies maintenance.
- Adjustable timer.

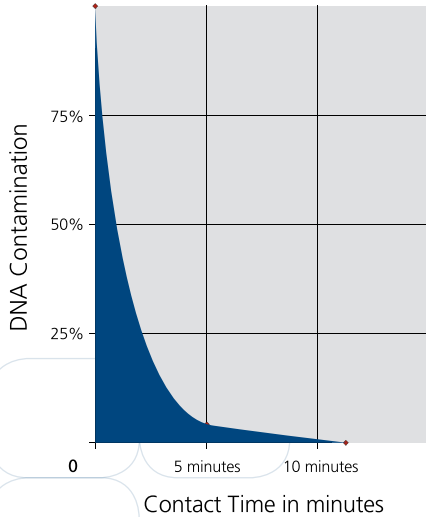


The Effect of UV on DNA Contamination

HEPA-Filtered Laminar Airflow

- Laminar, not turbulent flow provides superior sample protection.

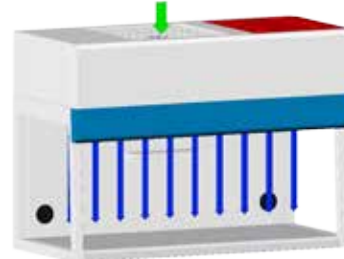

Laminar Flow (Esco PCR cabinets)

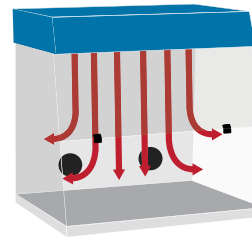

Turbulent Flow

Note: many competitors' cabinets have turbulent flow

- ISO Class 4 air cleanliness achieving clean air environment which is crucial to minimise the possibility of contaminating the precious samples - a key advantage against carrying out the critical procedure in open air or dead air box.
- Minipleat, separatorless HEPA filters, tested to a typical efficiency of $\geq 99.99\%$ for 0.3 micron particles are superior to conventional aluminium separator HEPA filters. Minipleat filters have a greater surface area and a longer service life, which reduce operating costs.

3

More Benefits

Easy-To-Use

- Timer is easy to adjust
- UV hour meter monitors bulb life
- Automatic decontamination for 0.9 m (3') and 1.2 m (4') models
 - Close sash: UV turns on automatically for decontamination
 - Open sash: airflow activates automatically

Safety

- UV interlock prevents UV exposure
- UL recognized electrical components
- UV filtering sash and side walls

Two-In-One


Can be used as a regular laminar flow cabinet and UV can also be used for decontamination of regular lab items.

Ergonomics

- Low noise
- Angled front to improve reaches
- Glass sides to prevent "boxed in" feeling
- Curved work surface front edge
- Powder-coated rear wall eliminates reflections
- Vertical clean airflow minimizes direct airflow towards operator, causing dry eyes



Energy Efficiency

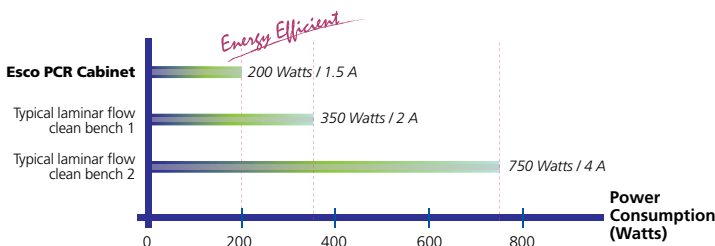

- Even more energy efficient than regular laminar flow clean benches
- UV timer shuts off UV after pre-set duration
- The backward curved wheel with external rotor motor delivers class-leading energy efficiency for lower operating costs

ESCO

WORLD CLASS. WORLDWIDE.

Esco PCR Cabinets

Provide Product Protection

Pre-Filters


An additional disposable pre-filter which is easily replaceable traps large particles in the inflow air prior to reaching the main filter, protecting against damage and prolonging filter life.

High Performance Fan System

German made ebm-papst® permanently lubricated, centrifugal motor/blowers with external rotor designs.


Motors selected for energy efficiency, compact design, and flat profile. Completely integrated assembly optimizes motor cooling.

All rotating parts balanced for smooth, quiet, vibration-free operation.

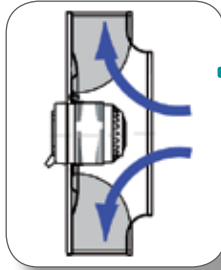

User Interface

An angled front, rounded work surface front edge, and glass sides promote ergonomics. The powder-coated work zone rear wall eliminates harsh reflections which may be associated with conventional stainless steel interiors. The vertical airflow design minimizes direct airflow which may lead to dry eyes and fatigue.

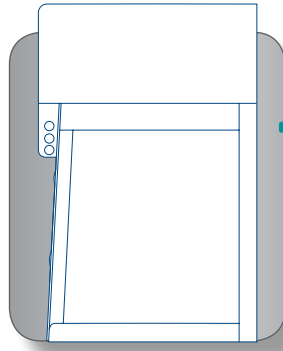

UV Decontamination System

Each cabinet includes a powerful, built-in, 253.7 nanometer, UV lamp to enable the work zone and materials to be decontaminated between experimental runs, thus preventing cross contamination.

The decontamination cycle time may be set with the UV timer, thus extending UV bulb life.



UV-filtering polycarbonate front and glass sides shield the user from accidental UV exposure. A proximity sensor ensures the frontal shield is lowered before allowing the UV to activate.

Air Cleanliness Standards

(ISO 14644-1, Air Cleanliness Particle Limits)
(No. of Particles / m³)

Particle Size (µm)	Cleanliness Class					
	Class 1	Class 2	Class 3	Class 4	Class 5	Class 6
0.1	10	100	1000	10000	100000	1000000
0.2	2	24	237	2370	23700	237000
0.3	-	10	102	1020	10200	102000
0.5	-	4	35	352	35200	35200
1.0	-	-	8	83	835	8320
5.0	-	-	-	-	29	293

Superior Air Cleanliness

Esco PCR cabinets provide ISO Class 4 air cleanliness within the work zone as per ISO 14644-1, significantly cleaner than the usual Class 5 classification on clean benches offered by the competition. The cabinets with the clean airflow are far better options than non ventilated box.

Esco PCR Cabinet,
Model PCR-4A_
with optional support stand



Control System




0.6 m (2') models are equipped with rocker switches for blower, light and UV

or




0.9 m (3') and 1.2 m (4') models are equipped with a microprocessor control system and soft touch controls for blower, light and UV.

Both models have a UV timer function; however, the microprocessor has additional program functions including automatic activation of UV when front shield is lowered, UV countdown on LCD display, total UV run hours, and more.

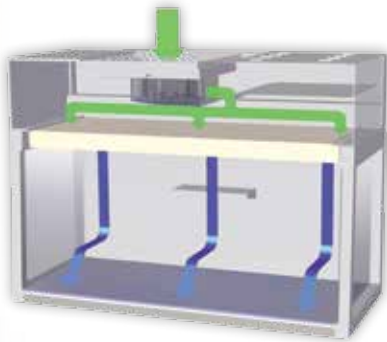

Proven Product Protection

Vertical laminar airflow with HEPA-filtration, $\geq 99.99\%$ at 0.3 microns, provides a sterile work space for PCR sample preparation.

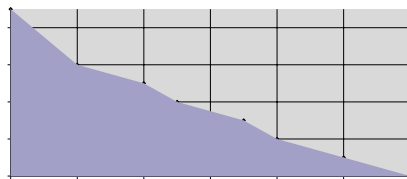

Built-In Protection

External surfaces are powder coated with Esco **ISOCIDE™** to eliminate 99.9% of surface bacteria within 24 hours of exposure.

Other Features

- All Esco products are manufactured for the most demanding clean air and containment requirements. All components are designed for maximum chemical resistance and enhanced durability for a long service life. The main body of the cabinet is constructed of industrial-grade electrogalvanized steel.
- One-piece formed stainless steel work surface with a rounded front edge is designed for maximum operator comfort.
- The lamp delivers uniform lighting to the work surface for greater comfort, reduced glare and improved productivity. The reliable lighting system is zero-flicker and instant start.
- Clean airflow achieves work area which is free from contaminants, crucial to ensure performance of the PCR process.
- Each PCR cabinet is individually factory tested for safety and performance in accordance with international standards.

Caution: PCR cabinets do not provide operator protection. They should not be used with applications involving unknown or hazardous agents.


General Specifications, Polymerase Chain Reaction Cabinets

Model	SCR-2A_	PCR-3A_	PCR-4A_				
Nominal Size	0.6 meters (2')	0.9 meters (3')	1.2 meters (4')				
External Dimensions (L x W x H)	730 x 617 x 950 mm 28.7" x 24.3" x 37.4"	1035 x 617 x 950 mm 40.7" x 24.3" x 37.4"	1340 x 617 x 950 mm 52.8" x 24.3" x 37.4"				
Internal Dimensions (L x W x H)	630 x 538 x 550 mm 24.8" x 21.2" x 21.7"	935 x 538 x 550 mm 36.8" x 21.2" x 21.7"	1240 x 538 x 550 mm 48.8" x 21.2" x 21.7"				
Laminar Airflow Velocity	0.275 - 0.325 m/s (54-64 fpm)	0.30 m/s (60 fpm)					
Pre-Filter	Washable polyurethane fibers with 85% arrestance						
Sound Emission*	<63 dBA	<56 dBA	<58 dBA				
Fluorescent Lamps Intensity	>800 Lux (>75 foot candles)	>975 Lux (>91 foot candles)	>1230 Lux (>114 foot candles)				
UV Lamp	253.7 nanometer 15-watt UV lamp		253.7 nanometer 30-watt UV lamp				
Controller	Rocker Switches	Esco Sentinel™ Microprocessor Control					
Construction	Main Body	Electrogalvanized steel with white oven-baked epoxy-polyester powder-coated finish. Coated with Esco Isocide™ antimicrobial coating					
	Work Zone	1.2mm (0.05") 18 gauge stainless steel Tyoe 304 with No.4 finish					
Electrical** (110-120V, AC, 60Hz, 1Ø)	Model	SCR-2A1	SCR-2A2	PCR-3A1	PCR-3A2	PCR-4A1	PCR-4A2
	Cabinet Full Load Amps (FLA)	1 A	1.5 A	0.75 A	1.1 A	1.8 A	3.8 A
	Cabinet Nominal Power	136 W	165 W	135 W	146 W	184 W	260 W
	Cabinet BTU****	464	563	461	498	628	887
Net Weight	76.5 kg (169 lbs)	99 kg (218 lbs)		115.4 kg (254 lbs)			
Shipping Weight***	100.5 kg (222 lbs)	123.3 kg (272 lbs)		139.8 kg (308 lbs)			
Shipping Dimensions, Maximum (L x W x H)***	850 x 730 x 1150 mm 33.5" x 28.7" x 45.3"	1130 x 730 x 1150 mm 44.5" x 28.7" x 45.3"		1420 x 730 x 1150 mm 55.9" x 28.7" x 45.3"			
Shipping Volume, Maximum***	0.71 m ³ (25 cu.ft.)	0.95 m ³ (34 cu.ft.)		1.19 m ³ (42 cu.ft.)			

6

- * Noise reading in open field condition / anechoic chamber
- ** Additional voltages may be available; contact Esco for ordering information.
- *** Cabinet only; excludes optional stand.
- **** Cabinet BTU = Cabinet nominal power x 3.41214.

Model PCR Cabinet Technical Specifications

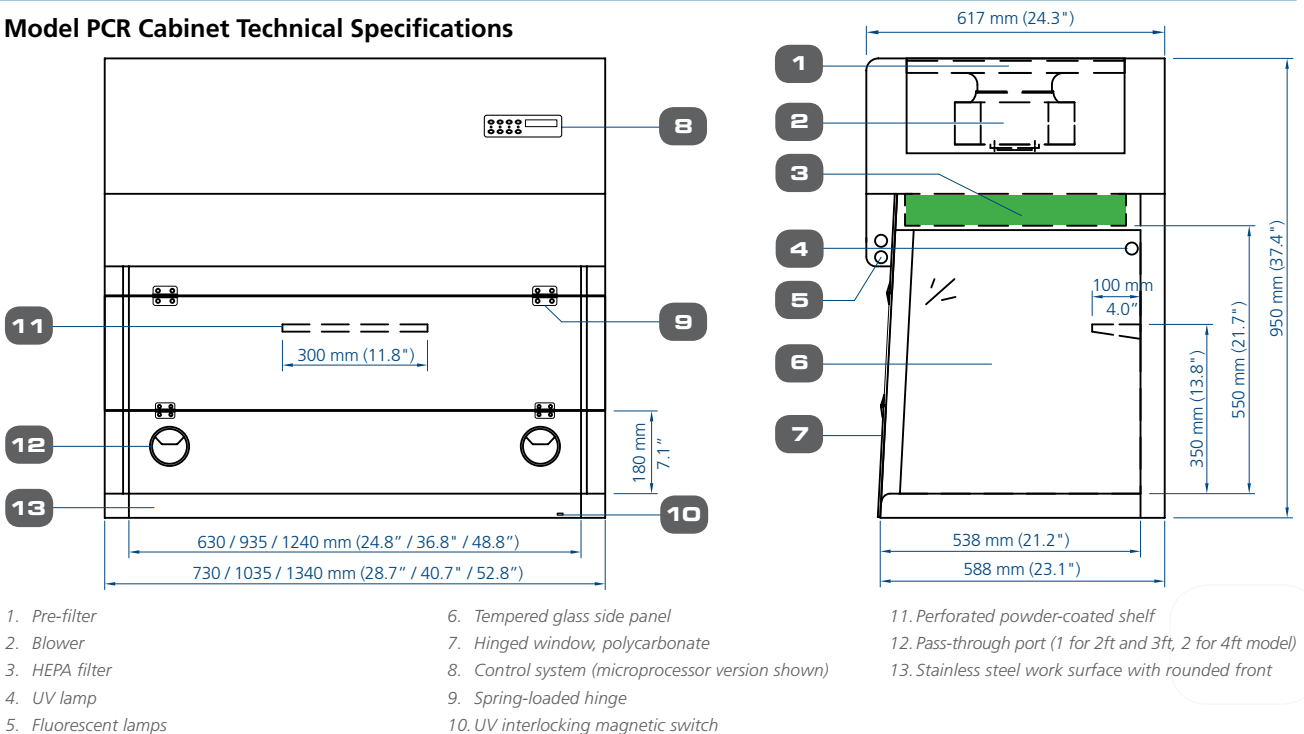

Standards Compliance	Cabinet Performance	Air Quality	Filtration	Electrical Safety
	IEST-RP-CC002, Worldwide	ISO 14644.1 Class 4, Worldwide AS 1386 Class 2.5, Australia JIS B9920 Class 4, Japan	IEST-RP-CC034, Worldwide IEST-RP-CC007, Worldwide IEST-RP-CC001, Worldwide EN1822 (H13), Europe	IEC 61010-1, Worldwide EN 61010-1, Europe UL-61010-1, USA CAN/ CSA 22.2 No. 61010-1

Accessories for PCR Cabinets



Support Stand with Caster Wheels (SPC)

- For 0.6 m (2'), 0.9 m (3') and 1.2 m (4') models
- Available in two standard heights: 711mm (28.0") or 860mm (34.0")
- Durable polyurethane caster wheels with 360 degree horizontal rotation
- Total brake system on front wheels
- Maximum weight supported: 600 kg (1323 lbs)



Support Stand with Leveling Feet (SAL)

- For 0.6 m (2'), 0.9 m (3') and 1.2 m (4') models
- Available in two standard sizes: 737 mm (29.0") or 864 mm (34.0"), ±38.1 mm (1.5")
- Maximum weight supported: 500 kg (1,100 lbs)



Telescoping Support Stands with Leveling Feet (STL)

- For 0.9 m (3') and 1.2 m (4') models
- Allow manual adjustment of the product height. The cabinet must be removed from a Telescoping Support Stand prior to adjustments
- Adjustable height range: 660-940 mm (26.0"-37.0"), adjustable in 25.4 mm (1.0") increments
- White oven-baked epoxy powder-coated finish
- Maximum weight supported: 600 kg (1323 lbs)



Telescoping Support Stands with Caster Wheels (STC)


- For 0.9 m (3') and 1.2 m (4') models
- Allow manual adjustment of the product height. The cabinet must be removed from a Telescoping Support Stand prior to adjustments
- Adjustable height range: 660-880 mm (26.0"-34.6"), adjustable in 25.4 mm (1.0") increments
- White oven-baked epoxy powder-coated finish
- total brake system on front wheels
- Maximum weight supported: 600 kg (1323 lbs)



Hydraulic Motorized Adjustable Support Stand with Casters (SPM)

- For 0.9 m (3') and 1.2 m (4') models
- Adjustable height range : 711-863 mm / (28.0"-34.0")
- Elevates to sitting or standing work surface height, Motorized electrically-adjustable
- Standard with caster wheels
- White oven-baked epoxy powder-coated finish
- Maximum weight supported: 500 kg (1100 lbs)

Note: Increases exterior dimensions

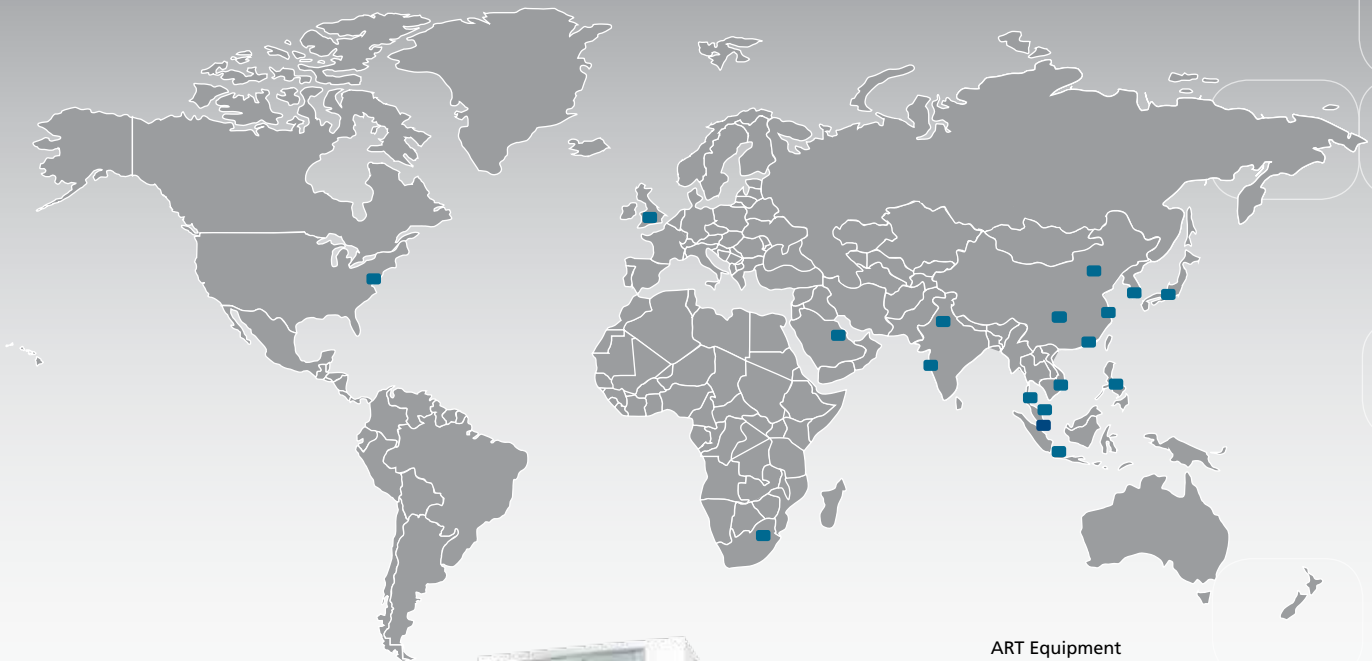

- ART Equipment
- Biological Safety Cabinets
- CO₂ Incubators
- Compounding Pharmacy Equipment
- Containment / Pharma Products
- Ductless Fume Hoods
- Freeze Dryer
- Lab Animal Research Products
- Laboratory Fume Hoods
- Laboratory Ovens and Incubators
- Laminar Flow Clean Benches
- PCR Cabinets
- PCR Thermal Cyclers
- Powder Weighing Balance Enclosures
- Ultra-low Freezers

The Esco Group of Companies is a global life sciences tools provider with sales in over 100 countries. The group is active in lab equipment, pharma equipment and medical devices. Manufacturing facilities are located in Asia and Europe. R&D is conducted worldwide spanning the US, Europe and Asia. Sales, service and marketing subsidiaries are located in 12 major markets including the US, UK, Singapore, Japan, China and India. Regional distribution centers are located in the US, UK, and Singapore.

Life Science • Chemical Research • Assisted Reproductive Technology (ART) • Pharmaceutical Equipment • General Equipment

ESCO

WORLD CLASS. WORLDWIDE.

Esco Technologies, Inc. • 2940 Turnpike Drive, Units 15-16 • Hatboro, PA 19040, USA
 Toll-Free USA and Canada 877-479-3726 • Tel 215-441-9661 • Fax 215-441-9660
us.escoglobal.com • usa@escoglobal.com

Esco Micro Pte. Ltd. • 21 Changi South Street 1 • Singapore 486 777
 Tel +65 6542 0833 • Fax +65 6542 6920 • mail@escoglobal.com
www.escoglobal.com

Esco Global Offices | Beijing, China | Chengdu, China | Kuala Lumpur, Malaysia | Manama, Bahrain | Guangzhou, China | Hanoi, Vietnam | Melaka, Malaysia | Mumbai, India | Philadelphia, PA, USA | Salisbury, UK | Shanghai, China | Seoul, Korea | Delhi, India | Osaka, Japan | Manila, Philippines | Midrand, South Africa | Jakarta, Indonesia | Singapore



9010022_PCR Cabinet_SCB & PCR_Combined Brochure_A4_9B3_0413
 Esco can accept no responsibility for possible errors in catalogues, brochures and other printed materials. Esco reserves the right to alter its products and specifications without notice. All trademarks and logos in this material are the property of Esco and the respective companies.